

Triclosan

• Intended Use

For detection of Triclosan and Triclosan methyl. Please refer to the attached specific procedures for water (groundwater, surface water, well water, effluent), and soil. Application procedures for other sample matrices can be obtained from Abraxis.

• Principle

The Abraxis Triclosan Assay applies the principles of enzyme linked immunosorbent assay (ELISA) to the determination of Triclosan and Triclosan methyl. The sample to be tested is added, along with paramagnetic particles attached with antibodies specific to Triclosan, to a disposable glass test tube and incubated for thirty (30) minutes. This is followed by the addition of a Triclosan enzyme conjugate. Both the Triclosan (which may be in the sample) and the enzyme labeled Triclosan derivative (the enzyme conjugate) compete for antibody binding sites on the magnetic particles. At the end of a thirty minute (30) incubation period, a magnetic field is applied to hold the paramagnetic particles (with Triclosan and labeled Triclosan analog bound to the antibodies on the particles, in proportion to their original concentration) in the tube and allow the unbound reagents to be decanted. After decanting, the particles are washed with Washing Solution.

The presence of Triclosan is detected by adding the enzyme substrate (hydrogen peroxide) and the chromogen (3,3',5,5'-tetramethylbenzidine). The enzyme-labeled Triclosan analog bound to the Triclosan antibody catalyzes the conversion of the substrate/ chromogen mixture to a colored product. After an incubation period of twenty (20) minutes, the reaction is stopped and stabilized by the addition of acid. Since the labeled Triclosan (conjugate) was in competition with the unlabeled Triclosan (sample) for the antibody sites, **the color developed is inversely proportional to the concentration of Triclosan in the sample.**

• Reagents

1. Triclosan Antibody Coupled Paramagnetic Particles

The Triclosan antibody (rabbit anti-Triclosan) is covalently bound to paramagnetic particles, which are suspended in buffered saline containing preservative and stabilizers.

100 test kit: one 60 mL vial

2. Triclosan Enzyme Conjugate

The horseradish peroxidase (HRP) labeled Triclosan analog is diluted in buffered saline containing preservative and stabilizers.

100 test kit: one 30 mL vial

3. Triclosan Standards

Five concentrations (0.025, 0.10, 1.0 ppb) of Triclosan in a methanolic/water solution with preservative and stabilizers. Each vial contains 2.0 mL.

4. Control

A concentration (approximately 0.50 ppb) of Triclosan in a methanolic/water solution containing preservative and stabilizers.

A 2.0 mL volume is supplied in one vial.

5. Diluent/Zero Standard

A methanolic/water solution containing preservative and stabilizers without any detectable Triclosan.

100 test kit: one 35 mL vial

6. Color Solution

A solution of hydrogen peroxide and 3,3',5,5'-tetramethylbenzidine in an organic base.

100 test kit: one 65 mL vial

7. Stopping Solution

A solution of diluted sulfuric acid (0.5%).

100 test kit: one 60 mL vial

8. Washing Solution

Preserved deionized water.

100 test kit: one 250 mL vial

9. Test Tubes

Glass tubes (36) are packed in a box.

100 test kit: three 36 tube boxes

• Reagent Storage and Stability

Store all reagents at 2-8°C. Do not freeze. Reagents may be used until the expiration date on the box. *The test tubes require no special storage condition and may be stored separately from the reagents to conserve refrigerator space.*

Consult state, local and federal regulations for proper disposal of all reagents.

• Materials Required but Not Provided

In addition to the reagents provided, the following items are essential for the performance of the test:

Pipets* Precision pipets capable of delivering 250 and 500 μ L and a 1.0 mL repeating pipet.

Vortex Mixer* Thermolyne Maxi Mix, Scientific Industries Vortex Genie, or equivalent

Magnetic Separation Rack*

Photometric Analyzer* capable of readings at 450 nm

* These items are available from Abraxis.

• Sample Information

Refer to sample preparation information contained under individual procedure (i.e. water, soil) or application notes.

Samples which have been preserved with monochloroacetic acid or other acids, should be neutralized with strong base e.g. 6N NaOH, prior to assay.

If the Triclosan concentration of a sample exceeds 1.0 ppb, the sample is subject to repeat testing using a diluted sample. A ten-fold or greater dilution of the sample is recommended with an appropriate amount of Diluent/Zero Standard or Sample Diluent. For example, in a separate test tube make a ten-fold dilution by adding 100 μ L of the sample to 900 μ L of Diluent/Zero Standard. Mix thoroughly before assaying. Perform the assay according to the Assay Procedure and obtain final results by multiplying the value obtained by the dilution factor e.g. 10.

The presence of the following substances up to 1000 ppm were found to have no significant effect on the Triclosan Assay results: copper, zinc, manganese, calcium, magnesium, sodium, phosphate, sulfate, FeSO₄, and nitrate. Humic acid and thiosulfate up to 100 ppm was found to have no significant effect.

• Reagent Preparation

All reagents must be allowed to come to room temperature and the antibody coupled paramagnetic particles should be mixed thoroughly before use.

• Procedural Notes and Precautions

As with all immunoassays, a consistent technique is the key to optimal performance. To obtain the greatest precision, be sure to treat each tube in an identical manner.

Add reagents directly to the bottom of the tube while **avoiding contact between the reagents and the pipet tip**. This will help assure consistent quantities of reagent in the test mixture.

Avoid cross-contaminations and carryover of reagents by using clean pipets for each sample addition and by avoiding contact between reagent droplets on the tubes and pipet tips.

Avoid foam formation during vortexing.

The magnetic separation rack consists of two parts: an upper rack which will securely hold the test tubes and a lower separator which contains the magnets used to attract the

antibody coupled paramagnetic particles. During incubations the upper rack is removed from the lower separator so that the paramagnetic particles remain suspended during the incubation. **For separation steps, the rack and the separator are combined to pull the paramagnetic particles to the sides of the tubes.**

To obtain optimum assay precision, it is important to perform the separation steps carefully and consistently. Decant the rack by slowly inverting away from the operator using a smooth turning action so the liquid flows consistently along only one side of the test tube. While still inverted, place the rack on an absorbent pad and allow to drain. Lifting the rack and replacing gently onto the pad several times will ensure complete removal of the liquid from the rim of the tube.

Mix the antibody coupled paramagnetic particles just prior to pipetting.

Standard and Control vials should remain capped when not in use, to prevent evaporation.

Do not use any reagents beyond their stated shelf life.

Avoid contact of Stopping Solution (sulfuric acid) with skin and mucous membranes. If this reagent comes in contact with skin, wash with water.

• Limitations

The Abraxis Triclosan Assay will detect Triclosan and Triclosan methyl to different degrees (refer to specificity table for data). The Abraxis Triclosan Assay kit provides screening results. As with any analytical technique (GC, HPLC, etc...) positive results requiring some action should be confirmed by an alternative method.

The total time required for pipetting the magnetic particles should be kept to two (2) minutes or less, therefore the total number of tubes that can be assayed in a run should be adjusted accordingly.

• Quality Control

A control solution at approximately 0.5 ppb of Triclosan is provided with the Abraxis Triclosan Assay kit. It is recommended that it be included in every run and treated in the same manner as unknown samples. Acceptable limits should be established by each laboratory.

• Assay Procedure

Read Reagent Preparation, Procedural Notes and Precautions before proceeding.

1. Perform the appropriate sample preparation according to the attached water or soil procedure. For any other sample matrices refer to specific procedures available from Abraxis.
2. Label **glass** test tubes for standards, control, and samples.

Tube Number	Contents of Tube
1,2	Diluent/Zero Standard, 0 ppb
3,4	Standard 1, 0.025 ppb
5, 6	Standard 3, 0.1 ppb
7, 8	Standard 4, 1.0 ppb
9	Control
10	Sample 1
11	Sample 2
12	Sample 3

3. Add 250 μ L of the appropriate standard, control, or sample.
4. Mix the Triclosan Antibody Coupled Paramagnetic Particles thoroughly and add 500 μ L to each tube.

5. Vortex for 1 to 2 seconds minimizing foaming.
6. Incubate for 30 minutes at room temperature.
7. Add 250 uL of Triclosan Enzyme Conjugate to each tube.
8. Vortex for 1 to 2 seconds minimizing foaming.
9. Incubate for 30 minutes at room temperature.
10. Separate in the Magnetic Separation Rack for **two (2) minutes**.
11. Decant and **gently** blot all tubes briefly in a consistent manner.
12. Add 1 mL of Washing Solution to each tube and **vortex tubes for 1-2 seconds**. Return tubes and allow to remain in the magnetic separation unit for **two (2) minutes**.
13. Decant and **gently** blot all tubes briefly in a consistent manner.
14. Repeat Steps 12 and 13 an additional time.
15. Remove the rack from the separator and add 500 uL of Color Solution to each tube.
16. Vortex for 1 to 2 seconds minimizing foaming.
17. Incubate for 20 minutes at room temperature.
18. Add 500 uL of Stopping Solution to each tube.
19. Add 1 mL Washing Solution to a clean test tube. Use as blank in Step 20.
20. Read results at 450 nm within 15 minutes after adding the Stopping Solution.

● Results

Manual Calculations

1. Calculate the mean absorbance value for each of the standards.
2. Calculate the %B/Bo for each standard by dividing the mean absorbance value for the standard by the mean absorbance value for the Diluent/Zero Standard.
3. Construct a standard curve by plotting the %B/Bo for each standard on vertical Ln (Y) axis versus the corresponding Triclosan concentration on horizontal Linear (X) axis on the graph paper provided.
4. %B/Bo for controls and samples will then yield levels in ppt of Triclosan by interpolation using the standard curve.
(Contact Abraxis for detailed application information on specific photometers.)

Photometric Analyzer

Some instrument manufacturers make available photometers allowing for calibration curves to be automatically calculated and stored. Refer to the instrument operating manual for detailed instructions. To obtain results from the Abraxis TRICLOSAN Assay on instruments allowing data transformation, the following parameter settings are recommended:

Data Reduct : Lin. Regression
 Xformation : Ln/Linear
 Read Mode : Absorbance
 Wavelength : 450 nm
 Units : PPT
 # Rgt Blk : 0

Calibrators:
 # of Cals : 4
 # of Repts : 2

Concentrations:
 #1: 0.00 PPT
 #2: 25 PPT
 #3: 100 PPT
 #4: 1000 PPT

Range : 20 - 1000
 Correlation : 0.990
 Rep. %CV : 10%

● Expected Results

Refer to the expected result section in the appropriate application note or procedure.

● Performance Data

Sensitivity

The Abraxis Triclosan Assay has an estimated minimum detectable concentration, based on a 90% B/Bo of 20 ppt (0.020 ppb). Refer to appropriate application notes or procedures for sensitivity in specific sample matrices.

Specificity

The cross-reactivity of the Abraxis Triclosan Assay for various related and unrelated compounds can be expressed as the least detectable dose (LDD) which is estimated at 90% B/Bo, or as the dose required to displace 50% (50% B/Bo).

Compound	LDD (ppb)	50% B/Bo (ppb)
Triclosan	0.020	0.250
Triclosan methyl	0.015	0.080
PBDE Congener 28	0.034	0.61
PBDE Congener 47	0.020	0.390
PBDE Congener 49	5.2	17.8
PBDE Congener 99	2.15	15.0
4'-OH-BDE-47	0.13	7.8
5-OH-BDE-47	0.15	5.6
6-OH-BDE-47	0.66	10.2
2,4,5-Tribromobiphenyl	> 100	> 100
2,4',5-Tribromobiphenyl	54	9,100
2,3,7,8-Tetrachloro-dibenzo-p-dioxin	> 100	> 100
T3	0.94	40
L-Thyroxine (T4)	340	700

The following compounds demonstrated no reactivity in the Triclosan Assay at concentrations up to 1,000 ppb: Biphenyl, 2,4-D.

● Assistance

For ordering or technical assistance contact:
 Abraxis LLC
 Sales Department
 54 Steamwhistle Drive
 Warminster, PA 18974
 (215) 357-3911 * Fax (215) 357-5232

● Ordering Information

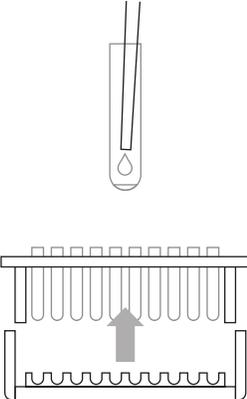
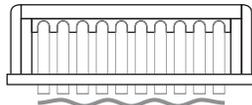
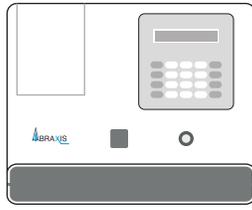
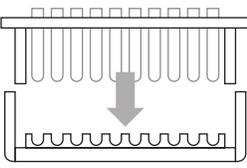
Abraxis Triclosan Assay Kit, 100T PN 530111
 Triclosan Sample Diluent PN 530112

● General Limited Warranty

Abraxis LLC warrants the products manufactured by the Company, against defects and workmanship when used in accordance with the applicable instructions for a period not to extend beyond the product's printed expiration date. **Abraxis makes no other warranty, expressed or implied. There is no warranty of merchantability or fitness for a particular purpose.**

022306

TRICLOSAN DETAILED FLOWCHART

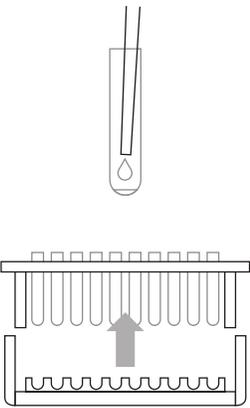
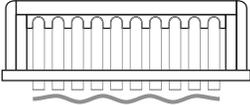
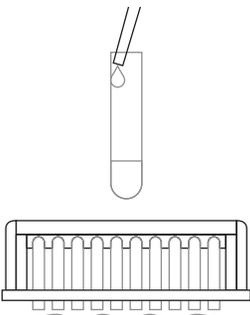
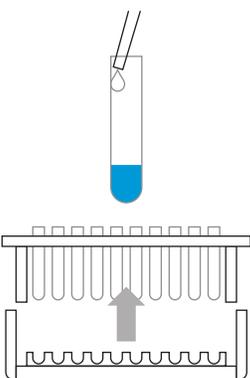
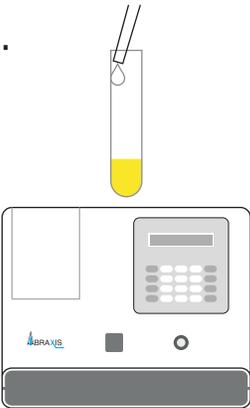
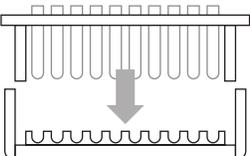
<p>1.</p>  <p><i>Remove</i> upper rack from magnetic base. Label test tubes for Standards, Control, and Samples.</p> <table border="1"> <thead> <tr> <th>Tube #</th> <th>Content</th> </tr> </thead> <tbody> <tr> <td>1, 2</td> <td>Diluent/Zero Standard 0 ppb</td> </tr> <tr> <td>3, 4</td> <td>Standard 1, 0.025 ppb</td> </tr> <tr> <td>5, 6</td> <td>Standard 2, 0.1 ppb</td> </tr> <tr> <td>7, 8</td> <td>Standard 3, 1.0 ppb</td> </tr> <tr> <td>9</td> <td>Control</td> </tr> <tr> <td>10</td> <td>Sample 1</td> </tr> <tr> <td>11</td> <td>Sample 2</td> </tr> <tr> <td>12</td> <td>Sample 3</td> </tr> </tbody> </table> <p><i>Add</i> 250 μL of either Standards, Control or Samples to the bottom of each test tube by inserting the pipette tip all the way into the bottom of the tube without touching the sides of the tube.</p>	Tube #	Content	1, 2	Diluent/Zero Standard 0 ppb	3, 4	Standard 1, 0.025 ppb	5, 6	Standard 2, 0.1 ppb	7, 8	Standard 3, 1.0 ppb	9	Control	10	Sample 1	11	Sample 2	12	Sample 3	<p>7.</p>  <p>Do not separate upper rack from lower base. Using a smooth motion, <i>invert</i> the combined rack assembly over a sink and pour out the tube contents; keep inverted and gently blot the test tube rims on several layers of paper toweling.</p>
Tube #	Content																		
1, 2	Diluent/Zero Standard 0 ppb																		
3, 4	Standard 1, 0.025 ppb																		
5, 6	Standard 2, 0.1 ppb																		
7, 8	Standard 3, 1.0 ppb																		
9	Control																		
10	Sample 1																		
11	Sample 2																		
12	Sample 3																		
<p>2.</p>  <p><i>Mix and Add</i> 500 μL of the Triclosan Antibody Coupled Paramagnetic Particles down the inside wall of each tube by aiming the pipet tip 1/4" to 1/2" below the tube rim without touching the rim or tube wall with the pipet tip; deliver liquid gently. <i>Vortex</i> for 1 to 2 seconds (at low speed to minimize foaming).</p>	<p>8.</p>  <p><i>Add</i> 1 mL of Washing Solution down the inside wall of each tube by using the technique described in Box 2. <i>Wait 2 minutes</i>. Using a smooth motion, invert the combined rack assembly over a sink and pour out the tube contents: keep inverted and gently blot the test tube rims on several layers of paper toweling. Repeat this step.</p>																		
<p>3.</p>  <p><i>React</i> 30 minutes at room temperature (15 °- 30°C).</p>	<p>9.</p>  <p><i>Lift</i> the upper rack (with its tubes) off the magnetic base; <i>add</i> 500 μL of Color Reagent down the inside wall of each tube by using the technique described in Box 2. <i>Vortex</i> for 1 to 2 seconds (at low speed to minimize foaming).</p>																		
<p>4.</p>  <p><i>Add</i> 250 μL of Triclosan Enzyme Conjugate down the inside wall of each tube by using the technique described in Box 2. <i>Vortex</i> for 1 to 2 seconds (at low speed to minimize foaming).</p>	<p>10.</p>  <p><i>React</i> for 20 minutes at room Temperature (15°- 30° C). During this period, add 1 mL of Washing Solution into a clean tube for use as an instrument blank in Step 10.</p>																		
<p>5.</p>  <p><i>React</i> 30 minutes at room temperature (15 °- 30°C).</p>	<p>11.</p>  <p><i>Add</i> 500 μL of Stopping Solution down the inside wall of each tube by using the technique previously Described. <i>Read</i> results at 450 nm within 15 minutes after adding the Stopping Solution. <i>Multiply</i> results of samples by the appropriate dilution factor (if any).</p> <p>[Safety Caution: Stopping Solution contains diluted sulfuric acid.]</p> 																		
<p>6.</p>  <p><i>Combine</i> the upper rack with the magnetic base; press all tubes into base; allow 2 minutes for the particles to separate.</p>																			

For Ordering or Technical Assistance Contact:
 ABRAXIS, LLC 54 Steamwhistle Drive, Warminster, PA 18974
 Phone: 215-357-3911 Fax: 215-357-5232
 Web: www.abraxiskits.com

Triclosan Magnetic Particle Kit Part # 530111, 100 Test



TRICLOSAN CONCISE FLOWCHART

<p>1.</p>  <p>Separate the rack.</p> <p>Add 250 μL of either Standards, Control or Samples to the bottom of each tube.</p>	<p>7.</p>  <p>Invert the combined rack.</p> <p>Blot gently.</p>
<p>2.</p>  <p>Mix the Triclosan Antibody coupled paramagnetic particles thoroughly and add 500 μL to each tube.</p> <p>Vortex for 1 to 2 seconds minimizing foaming.</p>	<p>8.</p>  <p>Add 1 mL of Washing Solution.</p> <p>Wait 2 minutes.</p> <p>Invert the combined rack.</p> <p>Blot gently.</p> <p>Repeat this step.</p>
<p>3.</p>  <p>Incubate for 30 minutes.</p>	<p>9.</p>  <p>Separate the rack.</p> <p>Add 500 μL of Color Reagent to each test tube.</p> <p>Vortex.</p>
<p>4.</p>  <p>Add 250 μL of mixed Triclosan Enzyme Conjugate to each test tube.</p>	<p>10.</p>  <p>Incubate for 20 minutes.</p> <p>Prepare blank.</p>
<p>5.</p>  <p>Incubate for 30 minutes.</p>	<p>11.</p>  <p>Add 500 μL of Stopping Solution to each test tube.</p> <p>Read OD 450</p>
<p>6.</p>  <p>Combine the rack and magnetic base.</p> <p>Seat all tubes.</p> <p>Wait 2 minutes.</p>	

For Ordering or Technical Assistance Contact:
 ABRAXIS, LLC 54 Steamwhistle Drive, Warminster, PA 18974
 Phone: 215-357-3911 Fax: 215-357-5232
 Web: www.abraxiskits.com

Triclosan Magnetic Particle Kit Part # 530111, 100 Test

